

# pH-Fix 0-14

Test strips for the determination of the pH-Value in the range of pH 0-14

## Contents:

100 Test strips  
1 Colour scale pH 0-14

## Instructions for use:

1. Dip test strip for 1 sec. with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with colour scale and read off the corresponding pH-value.

## Note:


If sample solution is weakly buffered leave test strip in solution until colour no longer changes.

## Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container below + 30 °C in a dry place.

## Disposal:

Please dispose used test strips in domestic waste.

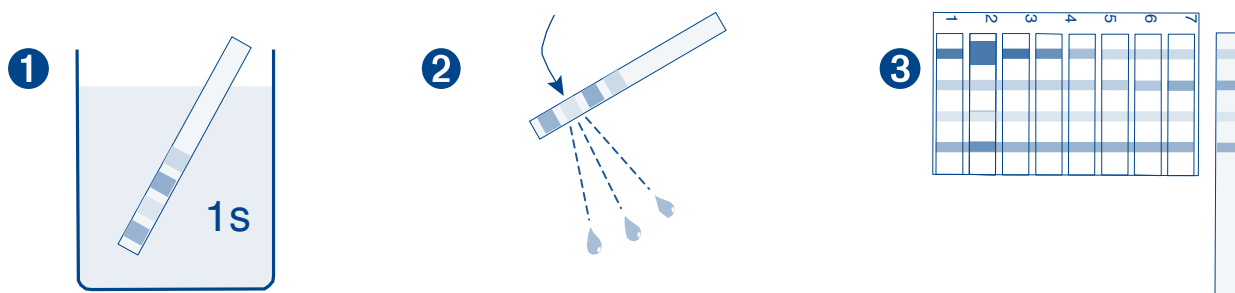

**MACHEREY-NAGEL GmbH & Co. KG** · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

Tel.: +49 24 21 969-0 · Fax: +49 24 21 969-199 · info@mn-net.com · [www.mn-net.com](http://www.mn-net.com)